

From Eye to Insight



Leica Microsystems CMS GmbH
Ernst-Leitz-Straße 17-37 | 35578 Wetzlar

Mateo FL RUO 1.2.2

Release Notes V 1.2.2.63

Pooja Dasgupta · Product Manager

pooja.dasgupta@leica-microsystems.com

Leica Microsystems CMS GmbH | Registered Office Wetzlar | Amtsgericht Wetzlar HRB 2432
Director: Dr. Annette Rinck | Gianmarco Cicuzza | Dr. Dirk Voelkel

Confidential - Company Proprietary

Mateo FL Release Notes

Abstract

This document describes the release of the Mateo FL software. You should read this document before installing your copy of this software.

All reasonable steps have been taken to ensure that this publication is correct and complete, but should any user be in doubt about any detail, clarification may be sought from Leica Microsystems (Suzhou) Technology Co., Ltd, or their accredited representative. The information in this document is subject to change without notice and should not be construed as a commitment by Leica Microsystems (Suzhou) Technology Co., Ltd. Leica Microsystems (Suzhou) Technology Co., Ltd accepts no responsibility for any errors that may appear in this document.

Copyright © 2026 Leica Microsystems (Suzhou) Technology Co., Ltd

All rights reserved. The contents of this publication may not be reproduced in any form, or communicated to a third party without prior written permission of Leica Microsystems (Suzhou) Technology Co., Ltd.

Due to a policy of continuous development, we reserve the right to change specifications without notice.

Date: July 2026

Printed in China

Introduction

This document describes the release of the Mateo FL software 1.2.2.63. Mateo FL is a fluorescence-equipped digital inverted microscope with precision automation tools and multi-channel fluorescence plus transmitted light imaging.

Microscope Support

Mateo FL

Mateo FL software

- When the new version is first run after installation or update, the EULA is displayed. The user reads and acknowledges the EULA terms before using the software.
- The user can stream live image, take pictures, record videos, control camera settings, repeat image settings, do measurements and annotations, check confluency, count cells, check transfection efficiency, save and share data to a USB, smart device or network folder.
- Supported languages: English and Chinese.

New features / changes:

This patch is to fix bugs and add improvements related to the CFR21p11 supporting features on Mateo FL.

- Admin can now be configured to prevent data collection and data generation activities.
- Dedicated user permission: System date and time, scale bar control, camera switch, image capture, contrast settings configurations are controlled through a dedicated user permission.
- Changes to system date and time, image format and filter cube settings are now captured in the Audit Trail, providing enhanced traceability of system configuration changes.
- Advanced Users can now be assigned Audit Trail Viewer rights.
- Configuration of network storage paths can be can now be enabled for Advanced Users.
- Admin-defined default settings are now enforced through role-based permissions. Settings such as Auto Save and Image Format can be locked to prevent unauthorized user modification, supporting consistency in data acquisition and storage workflows.

Software compatibility:

- The software runs only on Mateo FL.
- For updates, please check the product page at www.leica-microsystems.com